

# IP address change

Online version: [https://wiki.advacam.cz/wiki/IP\\_address\\_change](https://wiki.advacam.cz/wiki/IP_address_change)

# Contents



□  
Video of IP reset

- Connect the device to power.
- After the device init (R1 is **Orange**, changes to **Red** and turns off), attach the magnet to magnetic sensor. When the magnetic field is recognized, R1 turns **Blue**.
- When the change mode is enabled, R1 changes to one of the colors indicating the actual settings (listed below).
- Use magnet for select the preferred settings. There are 3 possibilities of the R1 color for different IP configurations:
  - **Red** = Auto IP
  - **Green** = DHCP
  - **Yellow** = Static IP (192.168.1.100)
- Wait for several seconds. R1 blinks several times and turns off. Now the setting is saved.

